

Whole Plasmid Sequencing – Express

Sample Preparation Guidelines

Service Overview

- Flat rate: \$4.99/sample
- Overnight data delivery
- Primer-free sequencing
- Low template input accepted (>10 µL)
- Texas customers only

Accepted Sample Types (Minimum Volume: 10 µL+)

- Plasmid DNA (5-100 ng/µL)
- RCA Product (10 ul)
- Agar Plate (No Agar plate accepted in Texas site)
- Bacterial Overnight Culture (10 ul)
- Colony Suspension (wash colony using 2 ul LB medium, transfer to 8x strips, add 20 ul medium, mix well, take 10 ul send to us; keep the rest volume as your backup)
- Glycerol Stock (10 ul)

Please Note, we do not keep nor return samples. Back up before sending out.

Sample Packing Instructions

Please use 8x strips. On tubes, only mark numbers according to order sheet is enough.

If shipping via public courier like UPS, Fedex, USPS etc:

- Place 8x strips in a strong enough container, eg. 50 ml tubes or card box (UPS overnight shipping offers free box). Unprotected tubes in envelop may results cap open, we experienced a lot. No ice necessary, try to avoid over heat by sunshine will be good enough.

Support

For any questions, please contact: jerry.li@quintarabio.com or text to 346-239-9971